



DBAP5

Manual version 1.16-EN
Firmware 1.5.2 or higher

5 Wavelengths Dual Beam Absorption Photometer



Via S. Andrea di Rovereto 33A CS - 16043 Chiavari (GE) Italy - Tel. +39 0185 318444

Email: len@len.it url: <http://www.len.it>



Summary

1. Scientific reasons	3
2. Instrument description	5
2.1. Pneumatic circuit	6
2.2. Measurement chamber	7
2.3. LED	8
2.4. Touch screen	8
2.5. Filter Tape	8
3. Measurement principle	9
3.1. Measurement cycle	11
3.2. Filter tape advancement	11
4. Local control (touch screen)	12
4.1. Measurement options	13
4.2. Calibrate	16
4.3. Open head	16
4.4. Download measure to USB stick	17
5. Remote control	18
5.1. Read Command	20
5.2. Write command (parameter setting)	21
5.3. Action command	22
5.4. Other commands	23
6. Changing the tape filter	23
APPENDICE A: data format	24
APPENDIX B: Flags	25
APPENDIX C: Electrical and mechanical specification	26
APPENDIX D: Normative	27
APPENDIX E: Poster PM2022	28

1. Scientific reasons

Atmospheric aerosol, both of natural and anthropogenic origin, regulates the earth global temperature by acting as a forcing on the radiative balance due to its optical properties, i.e., the ability to absorb and diffuse sunlight.

The absorption of solar radiation act as a regulator of the global terrestrial temperature, in fact the heating to the atmosphere and the cooling to the Earth surface caused by this absorption are hypothesized to have significant climate impacts¹, therefore reliable observations of the aerosol absorption coefficients (K_{ab}) are crucial to quantify the radiative forcing of the climate, and so aerosol absorption measurement are recommended for all station in the GAW network.

From the absorption coefficients it is possible to retrieve the equivalent black carbon (EBC) concentration, using the adequate value of Mass Absorption Coefficient (MAC) in **Eq. 1**

$$EBC = \frac{K_{ab}^{\lambda}}{MAC^{\lambda}}$$

Eq. 1

EBC can have effect not only on air quality but also on human health, as highlighted by the WMO report².

A key factor in better understanding of the aerosol composition is the spectral dependency of absorption coefficient given by the Angstrom Absorption Exponent (AAE) defined as **Eq. 2**

$$K_{ab}^{\lambda} \propto \lambda^{-AAE}$$

Eq. 2

The values of AAE between 870 nm and 420 nm, calculated from **Eq. 3** as

$$AAE_{420-870} = \left(\frac{\ln(K_{ab}^{870} / K_{ab}^{420})}{\ln(870/420)} \right)$$

Eq. 3

¹ *Particulate absorption of solar radiation: anthropogenic vs. dust. C. Wang et al 2009 Atmos. Chem. Phys.*

² *Health effects of black carbon 2012 World Health Organization.*

This is widely used to characterize the particle composition, as reported in³ where AERONET AOD dataset has been investigated. From this work, Figure shows that different AAE values are related to different aerosol composition.

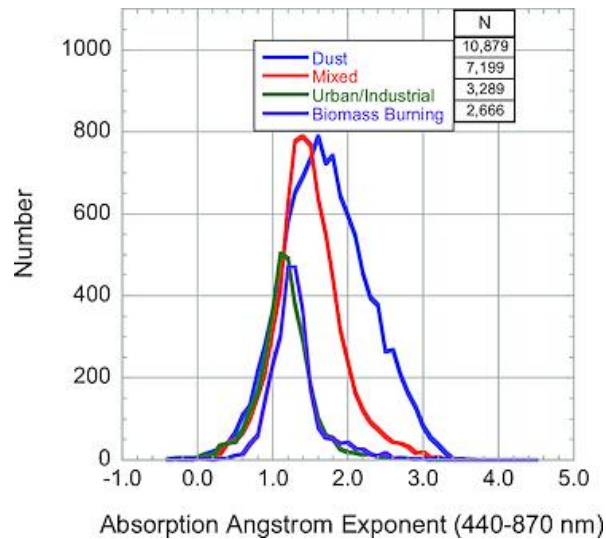


Figure 1

³ Analysis of AERONET aerosol absorption properties and classifications representative of aerosol source regions. D. M. Gilles et al. 2012 J. Geophys. Res.

2. Instrument description



Figure 2

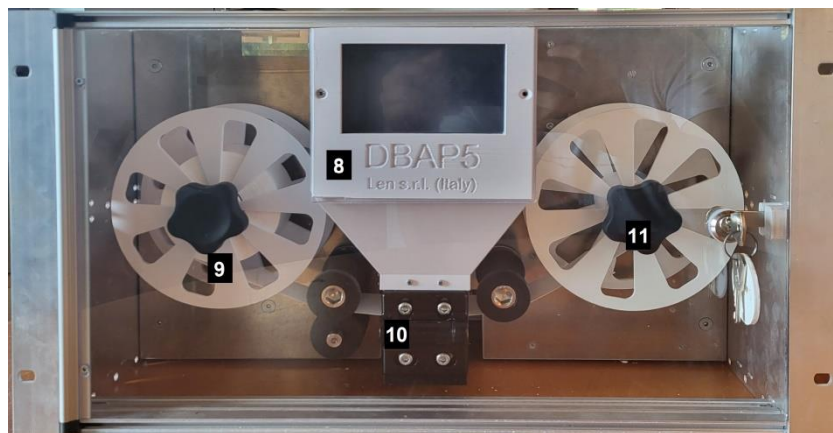


Figure 3

- 1 Air inlet
- 2 Air outlet
- 3 Power plug and switch
- 4 Ethernet plug RJ45
- 5 Female RS232 db9 port
- 6 USB port
- 7 Cooling fan
- 8 LCD screen with touch screen
- 9 Filter tape
- 10 Measurement chamber
- 11 Plexiglas cover lock

2.1. Pneumatic circuit

Figure 4 shows the pneumatic circuit of the DBAP5. Air flow entering the measuring chamber (1), after the measuring channel the flow is measured through the Venturi tube (2). The flow then enters the reference channel. Any possible flow fluctuation due to the pump (5) is attenuated by an expansion lung (4). The air flow is then evacuated through the instrument outlet.

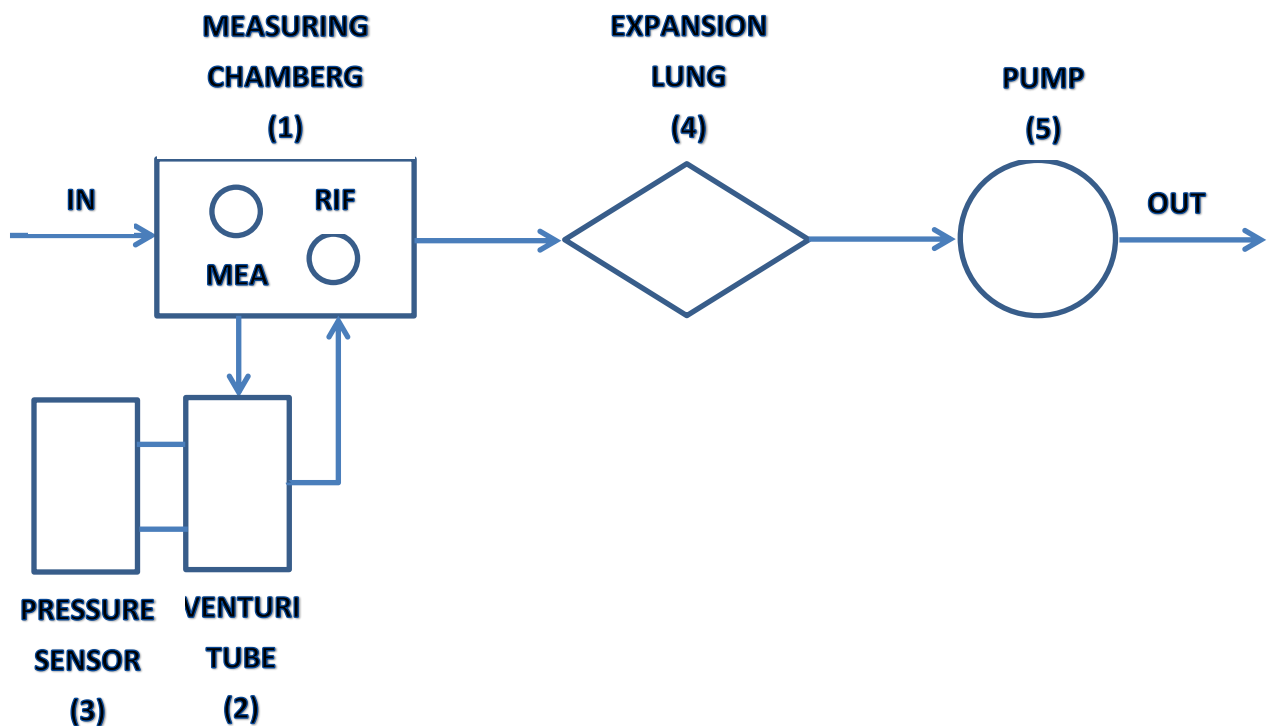


Figure 4

Air flow in the Venturi tube is measured using Eq. 4

$$Q = a + b\sqrt{dP * 100}$$

Eq. 4

In which Q is the air flow in l/min and dP is the Venturi tube's differential pressure. The factor 100 is necessary to convert dP from millibar (unit used by instrument) to Pascal.

Coefficients a and b mainly depend on the geometric characteristics of the tube.

In DBAP5 they are set following a factory calibration, however they can be changed following a further calibration performed by user⁴.

Sampled air relative humidity (RH) is measured between the inlet and the measurement chamber, while sampled air temperature is measured inside the Venturi tube. Both air RH and temperature offset can be adjusted by user. Atmospheric pressure is also recorded.

2.2. Measurement chamber

The measuring chamber (**Figure 5**), designed according to dual beam technology, has two distinct cells (reference and measurement) in which the light generated by the LED is conveyed using an optical fiber. This technique, ensuring a continuous comparison with the uncontaminated filter during the measurement, cuts any errors due to possible changes in the optical properties of the filter due, for example, to changes in humidity and temperature. The light intensity is in fact measured downstream of the filter in both cells, supplying the transmittance at each instant of time and for each wavelength, the variation of which over time supplies the absorption coefficient (paragraph 3).

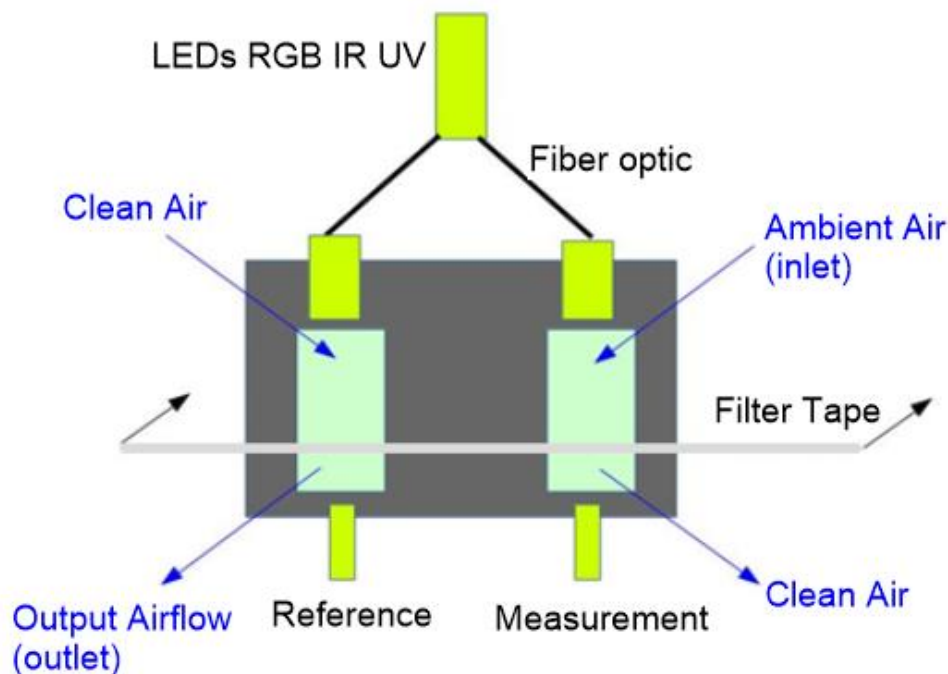


Figure 5

⁴ See paragraph 5.3 for parameters setting

Ambient air enters the upper part of the measuring cell, passes through the filter, depositing the particulate load on it, and exits the lower part of the cell.

After passing through the filter, the air is free of aerosols and enters the reference cell from above.

The developed measuring chamber ensures that the filter is subjected to the same pressures to minimize the possible difference in the optical path of the light inside the filter due to different deformations of the filter itself.

2.3. LED

Light sources are at the following wavelength:

- IR: 870 nm
- RED: 634 nm
- GREEN: 522 nm
- BLUE: 465 nm
- UV: 420 nm

2.4. Touch screen

The instrument display is a touch screen that allow the user to perform basic operation on the instrument and control the measured data. The use of touch screen is described in paragraph 4.

2.5. Filter Tape

The filter tape used for the measurement is made of fiberglass (GF 10 width 40 mm, retention degree 99.98%). To keep measure efficiency, the filter must be advanced when particles load became too intensive, thus the advancement is driven by transmittance value, as described in paragraph 3.2, that can be set by the user.

For the filter replacement procedure, refer to paragraph 6.

3. Measurement principle

The method for finding the absorption coefficient for each wavelength, and then the equivalent black carbon, is based on the measurement of the change in filter transmittance as particulate matter accumulates on it.

Using the 'dual beam' technology (paragraph 2) the transmittance (τ) is defined for each wavelength as the ratio between the light intensity measured in the measuring cell (I_M) and in the reference cell (I_R), transmittance which is normalized to 1 in the first measurement cycle ($t = 0$).

$$\tau^\lambda(t) = \frac{I_m^\lambda(t)/I_r^\lambda(t)}{I_m^\lambda(t=0)/I_r^\lambda(t=0)}$$

Eq. 5

the attenuation coefficient for each wavelength (K_{att}) in unit of m^{-1} is calculated from the filter transmittance variation over a time interval following **Eq. 6**

$$K_{att}^\lambda(t) = \frac{A}{Qdt} \ln \left(\frac{\tau^\lambda(t-dt)}{\tau^\lambda(t)} \right)$$

Eq. 6

In which A is the area of the filter where particles are collected (spot area) in m^2 , Q is the air flow passing through the filter in m^3/s and dt the time interval in seconds between the two measurements, which is set to 60 s.

The absorption coefficients are then calculated from the attenuation coefficients taking in account of two effects on the aerosol optical properties due to the filter media, using the correction factor **Eq. 7**

$$F(\tau) = \frac{1}{C(a + b\tau)}$$

Eq. 7

where the coefficient C depends mainly on the type of filter and considers the increase in absorption due to multiple scattering within the filter fiber.

F(t) depends linearly on the transmittance and describe the shadow effect that leads to an underestimation of absorption due to the growth of the particulate load on the filter itself.

In **Eq. 7** C, a and b can be set by the user, and by default they are set as:

- a = 0.531
- b = 0.610
- C = 2.3+0.0003λ[nm]

the absorption coefficients (K_{ab}^{λ}) are therefore calculated as:

$$K_{ab}^{\lambda} = K_{att}^{\lambda} F(\tau)$$

Eq. 8

The Equivalent Black Carbon (EBC) is derived from the absorption coefficients using **Eq. 1**.

DBAP5 gives the value of EBC calculated at 870 nm (IR), using a MAC of 6.17e-6, value that can be set by the user following the scientific upgrade. The DBAP5 also allows the user the possibility to set a MAC value for each wavelength to get EBC for all the 5 colours.

$$BC = \frac{K_{ab}^{\lambda=870}}{MAC^{\lambda=870}} [\mu g/m^3]$$

Eq. 9

To minimize measurement noise, especially in clean environment, the user can decide to receive the absorption coefficient as simple moving average (SMA) over a time (P) of a maximum of 10 minutes.

$$K_{ab}^{\lambda} = \sum_{i=-P}^{i=0} K_{ab}^{\lambda}(i)$$

Eq. 10

where i=0 represent the last received data.

3.1. Measurement cycle

After closing the measurement chamber, the pump is turned on and the dark current noise is measured with LED off. The background noise thus obtained will be subtracted from the next light intensity measurements. This phase has a total duration of 5 second, as 1 seconds are used for each color.

Subsequently all five transmittances are normalized to 1. This phase has a total duration of 60 seconds, as 12 seconds are used for normalization for each color.

After these first two phases, the real measurement cycle begins: transmittance (**Eq. 5**) is measured every second for each wavelength and integrated over a minute to perform the calculation of the absorption coefficient in m^{-1} (**Eq. 8**) and of the equivalent black carbon in $\mu\text{g}/\text{m}^3$ (**Eq. 9**).

After P minutes (chosen SMA interval) the first value of the absorption coefficients and EBC is supplied for each of the 5 wavelengths, values that are updated every minute.

3.2. Filter tape advancement

As the amount of absorbing matter collected on the filter reaches a critical value, the light intensity variation becomes very weak leading big uncertainty in the measure. Thus, the filter tape is advanced when the transmittance values reaches the value set in the configuration for each color, default values are 0.6 for UV, other colors set to 0.0 (disabled) .

When this transmittance threshold is reached the measure is stopped, the pump gets turned off and the measurement chamber lifts its upper part to let the filter tape pass inside. After the filter is in the new position the measurement chamber closes and the measure starts again.

The user can in any case change the default transmittance values. In the case of limits set to multiple colors, the first threshold reached will shift the filter.

4. Local control (touch screen)

After the instrument has been turned on, it is able to receive remote commands (RS22, telnet), and after touching the screen the instrument can receive local commands, following the menu shown in **Figure 6**. The last button (DOWNLOAD MEASURES TO USB STICK) appears only if a USB stick is inserted in the instrument.

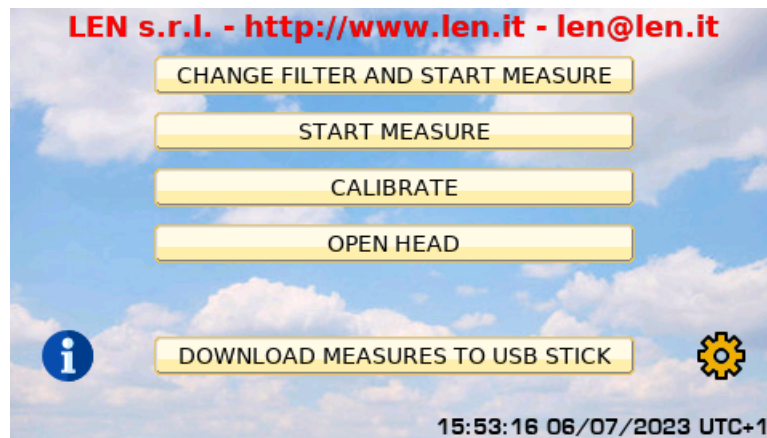


Figure 6

Choosing the maintenance icon on the bottom-right corner of the screen it is possible to change language (Italian, English or French) used on the touch screen.

The icon information shows a new screen with some of the configuration of the instrument, as shown in **Figure 7**

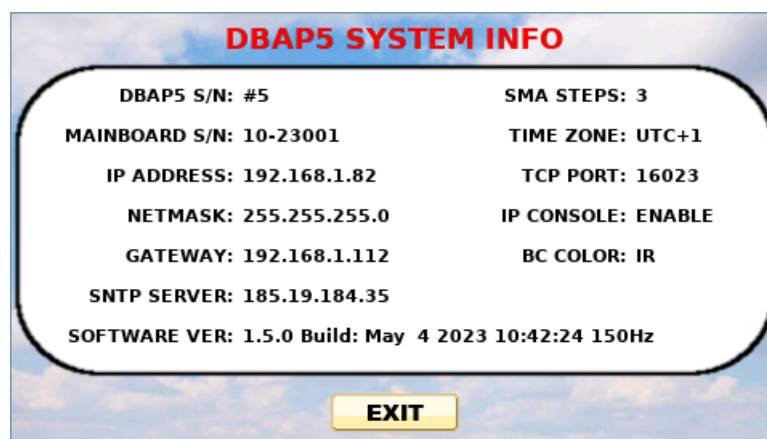


Figure 7

4.1. Measurement options

In the main menu (**Figure 6**) the top first two options (“CHANGE FILTER AND START MEASURES” and “START MEASURES”) activate the measures.

By choosing the first choice the filter tape will advance and the measurement cycle start, and by choosing the second choice the measurement cycle will start without changing the filter position.

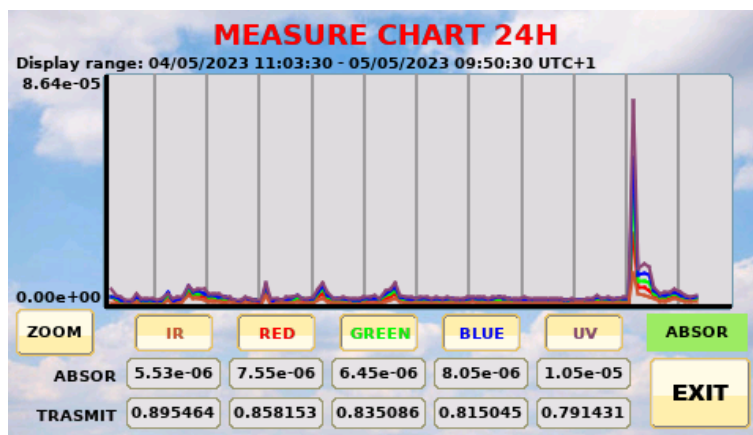


Figure 8

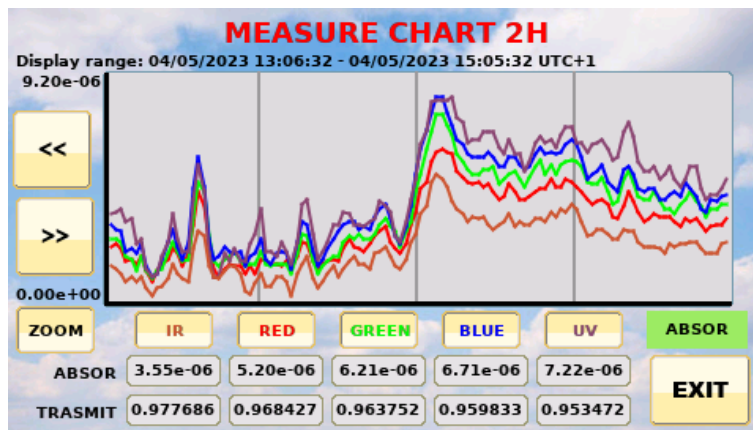


Figure 9

During measurement operations, the display will show the time series of the absorption coefficients measured for all wavelengths during the time interval showed at the top of the screen (Range display time). The vertical grid shows 2-hour intervals (or 30 minutes in case of activated zoom). The last measured values of absorption coefficient (in m^{-1}) and transmittance are reported, for each wavelength, in the boxes at the bottom of the screen. By clicking on the options that show the name of the colours, you can choose whether to display the absorption coefficient measured at the relative

wavelength. 1 to 5 curves can be displayed at the same time. However, at least one time series is always displayed.

The ZOOM button allows the last 2 hours of measurement to be displayed on the screen instead of the last 24 hours and vice versa. If zoom is enabled the [] navigation button displayed to the right of the chart ordinate moves the zoom time window one hour back and the [>>] button one hour forward.

The EXIT button can be used to stop the measurement.

The green bottom on the bottom right allows to switch between graphical visualization of absorption coefficients (ABSOR **Figure 8** and **Figure 9**), Angstrom Absorption Exponent (AAE **Figure 10**), transmittances (TRANSMIT **Figure 11**) and a screen reporting numerical values (VALUES **Figure 12**), equivalent black carbon (BC **Figure 13**).

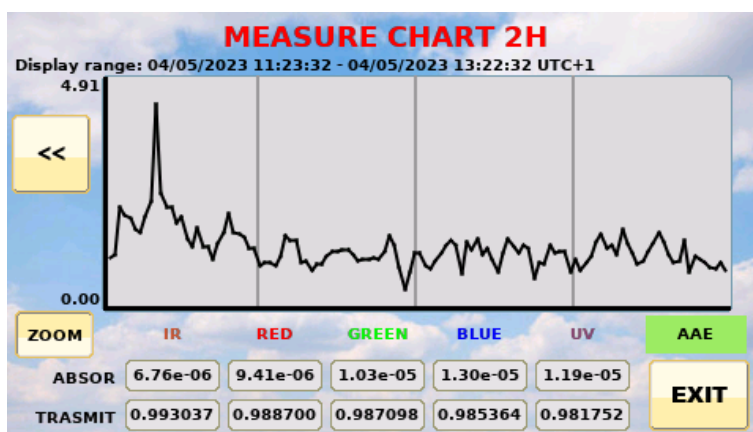


Figure 10

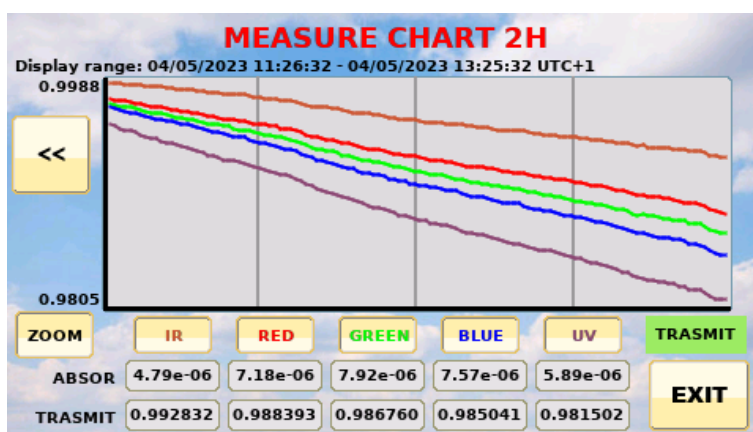


Figure 11

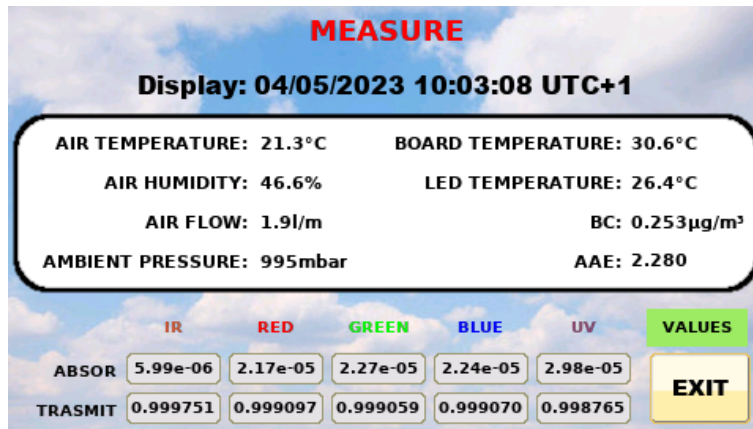


Figure 12

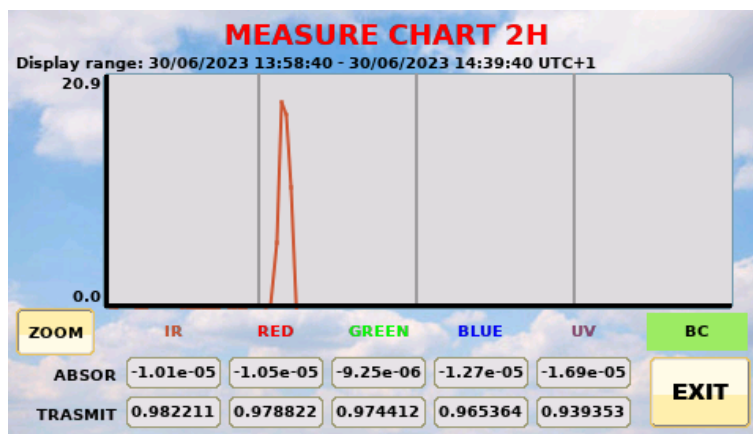


Figure 13

4.2. Calibrate

The CALIBRATE button put the instrument in the calibration mode, the mode that is mainly used by the manufacturer. In this mode, the screen will display the signal received by the two detectors of the measurement chamber (MEASURE and REFERENCE) together with the related transmittance to check the signal stability (**Figure 14**).

The screen shows the result of the signal in the selected wavelength, choosing NO LIGHT the LED will be turned off and the dark current measurement is performed.

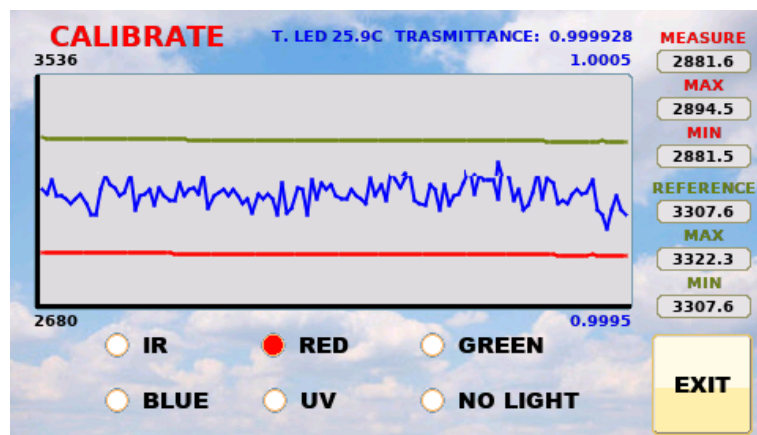


Figure 14

4.3. Open head

The OPEN HEAD button forces the measurement chamber to open, and the screen of **Figure 15** to appear on the display. From this new menu, it's possible to advance the filter (by choosing FORWARD REEL), increase the opening of the chamber (OPEN HEAD) or close the measurement chamber (CLOSE HEAD).



Figure 15

4.4. Download measure to USB stick

This choice appears only if a USB mass storage is plugged in the USB port of the DBAP and measure is not active. Choosing this option, a new screen (Figure 16) allows to decide the period of data that must discharged.

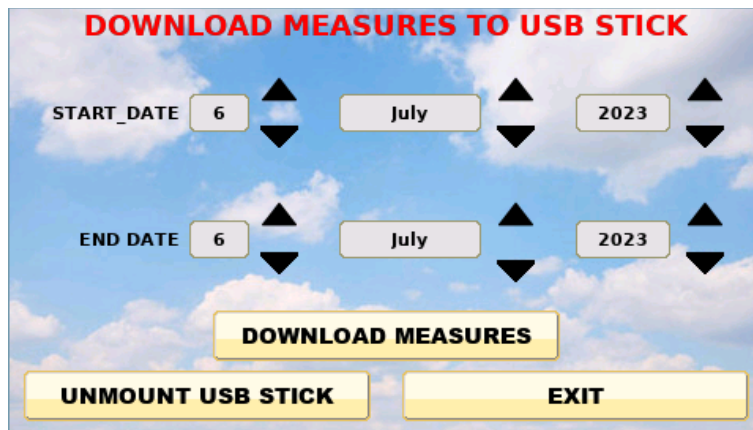


Figure 16

Note: to save data to USB measurement must stopped!

5. Remote control

Using the communication protocol via RS232 or telnet allows the user to gain greater control of the instrument. The RS232 port has the following parameters: 57600, 8, N, 1. The telnet uses port 16023 as the default port, but it can be changed with the CHANGE_PORT command.

Network settings: must be provided by the user exclusively via the serial link (command: CHANGE_IP and CHANGE_GATEWAY, paragraph 0).

To visualize on a computer terminal all possible commands type “?” .

Command	Description
?	Show Help Menu
ADMIN	Switch to admin menu. Reserved for the producer – need a password.
CHANGE_FILTER	Change filter
CHANGE_GATEWAY	Change IP gateway (CHANGE_GATEWAY IPSUB1.IPSUB2.IPSUB3.IPSUB4)
CHANGE_IP	Change IP address and netmask (CHANGE_IP IPSUB1.IPSUB2.IPSUB3.IPSUB4 NETSUB1.NETSUB2.NETSUB3.NETSUB4)
CHANGE_PORT	Change TCP remote console port (CHANGE_PORT NUM_PORT)
CHANGE_SNTP_SERVER	Change SNTP IP SERVER (CHANGE_SNTP_SERVER IPSUB1.IPSUB2.IPSUB3.IPSUB4). With no arguments disabled SNTP
CLOSE_HEAD	Close head
CLOSE_REMOTE_SESSION	Close remote session (console telnet)
DATE	Display date and time
DELETE_ALL_MEASURES	Delete all measures
DELETE_DATE_MEASURES	Delete all measures of that day (DELETE_DATE_MEASURES DD/MM/YYYY)
DELETE_LOGS	Delete all logs
DISABLE_FLUXMETER	Disable the fluxmeter The flow is fixed to user default
DISABLE_PUMP	Disable the pump in the measure. The flow is fixed to user default
DISABLE_REMOTE_CONSOLE	Disable remote console
ENABLE_FLUXMETER	Enable the fluxmeter.
ENABLE_PUMP	Enable the pump in the measure.
ENABLE_REMOTE_CONSOLE	Enable remote console
FORWARD_REEL	Forward reel
GET_DATE_MEASURES	Print measures of date (GET_DATE_MEASURES DD/MM/YYYY [hh:mm]).
GET_DATE_RANGE_MEASURES	Print measures date range (GET_DATE_MEASURES DD/MM/YYYY DD/MM/YYYY).
GET_HEADER_MEASURE	Print header measures.
GET_LAST_MEASURES	Print last measures (Max. 1 day). Without argument show only last
LED_BLUE	Switch on blue led
LED_GREEN	Switch on green led
LED_IR	Switch on IR led
LED_NONE	Switch off all led

Command	Description
LED_RED	Switch on red led
LED_UV	Switch on UV led
OPEN_HEAD	Open head
READ_AIR_HUMI	Read air humidity
READ_AIR_TEMP	Read flux temperature
READ_ENVIRO_PRESSURE	Read environmental pressure
READ_FLUX_L_MINUTE	Read flux l/min
READ_FLUX_PRESSURE	Read flux pressure
READ_INT_TEMP	Read internal temperature
READ_LED_TEMP	Read led temperature
SET_AIR_HUMI_OFFSET	Set air humidity offset (range -10 to +10)
SET_AIR_TEMP_OFFSET	Set air temperature offset (range -10 to +10)
SET_BC_CONSTANT	Set constant to calculate black carbon (syntax SET_BC_CONSTANT <COLOR> value)
SET_DATE	Set date and time (hh:mm:ss DD/MM/YYYY)
SET_DEFAULT_FLUX	Set default flux for no pump or no fluxmeter
SET_FACTORY_CONFIG	Set factory configuration and restart
SET_FILTER_CORRECTION_CONST	Set filter correction const (syntax SET_FILTER_CORRECTION_CONST <COLOR> <CONST_A> <CONST_B> <CONST_C>)
SET_FLOW_CONST_A	Set flow const A
SET_FLOW_CONST_B	Set flow const B
SET_LEVELS_FILTER_CHANGE	Transmittance level filter change for colors. (IR_TR RED_TR GREEN_TR BLUE_TR UV_TR)
SET_SMA_STEPS	Set Simple Moving Average steps. 1 to 10 measures
SET_SPOT_AREA	Set spot area
SET_TIME_ZONE	Set UTC time zone
SHOW_CONFIG	Show configuration parameters
SHOW_DATE_MEASURES	Show date measure present on sd (optional parameters DATE NUM_DATE_MEASURES).
SHOW_LOG_GEN	Show generic logs. (Optional argument last number of logs).
SHOW_LOG_SYS	Show system logs. (Optional argument last number of logs).
SHOW_STATUS	Show DBAP status
SHOW_VER	Show software version
SPD	Set pump duty cycle
START_CALIBRATION	Measure and reference every second update
START_MEASURE	Start measure. Option PUMP enable at once the pump
START_PUMP	Start pump
START_STREAMING	Start streaming measure
STOP_CALIBRATION	Stop calibration
STOP_MEASURE	Stop measure (switch off clock, led and pump)
STOP_PUMP	Stop pump
STOP_STREAMING	Stop streaming measure



Commands are not case sensitive, and they must be followed by the \n\r characters.

DBAP5 commands can be divided in Action, Write and Read commands. Action commands are used for activating some operation, Write commands for parameters setting and Read commands to receive information and data for the instrument.

5.1. Read Command

Command	
?	Show Help Menu
DATE	Show instrument date and time
GET_DATE_MEASURES	Download data of the given day starting from the given time (time is optional), data format is reported in appendix A.
GET_DATE_RANGE_MEASURES	Download all data in the given dates range (GET_DATE_MEASURES DD/MM/YYYY DD/MM/YYYY).
GET_HEADER_MEASURE	Download the header as reported in Appendix A.
GET_LAST_MEASURES	Show last measurement data, it can be given with an integer (n) less than 1440 as argument to get last n measurement data, in is missing only the last measure is downloaded. Data format is reported in appendix A.
READ_AIR_HUMI	Show the air sampled air temperature [%].
READ_AIR_TEMP	Show the temperature [Celsius degrees] inside the Venturi tube.
READ_ENVIRO_PRESSURE	Show enviroment pressure [mbar].
READ_FLUX_L_MINUTE	Show flow [l/min].
READ_FLUX_PRESSURE	Show the differential pressure [mbar] in the Venturi tube.
READ_INT_TEMP	Show the internal temperature [Celsius degrees].
READ_LED_TEMP	Show led temperature [Celsius degrees].
SHOW_CONFIG	Show the parameter setting on the instrument.
SHOW_DATE_MEASURES	Show the date of the measures saved in the instrument memory, if used with argument it will show the date starting from the given date.
SHOW_LOG_GEN	Shows the instruments generic logs (maximum last 1000 events).
SHOW_LOG_SYS	Shows the instruments system logs (maximum last 1000 events).
SHOW_STATUS	Shows instrument status and errors.
SHOW_VER	Shows the firmware and bootloader version.
START_STREAMING	Start data streaming over RS232 or telnet. Data format is reported in appendix A
STOP_STREAMING	Stop data streaming

5.2. Write command (parameter setting)

Command	Description
CHANGE_GATEWAY	Only over RS232 Set the IP gateway.
CHANGE_IP	Only over RS232 Set IP and netmask.
CHANGE_PORT	Set the TCP port for the communication over telnet.
CHANGE_SNTP_SERVER	Set the NTP server for time synchronization. Used without parameter will disable the time synchronization.
SET_AIR_HUMI_OFFSET	Set the offset for the air humidity sensor by adding the given value.
SET_AIR_TEMP_OFFSET	Set the offset for the air temperature sensor by adding the given value.
SET_BC_CONSTANT	Set the color and the related MAE of Eq. 8 to calculate the EBC.
SET_DATE	Set instrument date and time.
SET_DEFAULT_FLUX	Set the fictitious flux value for operations without active pump.
SET_FACTORY_CONFIG	Set instrument factory configuration (appendix B).
SET_FILTER_CORRECTION_CONST	Set the coefficient for the filter correction function Eq. 7.
SET_FLOW_CONST_A	Set coefficients a for the air flow calibration Eq. 4 .
SET_FLOW_CONST_B	Set coefficients b for the air flow calibration Eq. 4 .
SET_LEVELS_FILTER_CHANGE	Set the transmittance threshold for advancing the filter tape (paragraph 3.2). Setting the value to 0 will disable the tape movement for the related colour.
SET_SMA_STEPS	Set the number of measures to use for the SMA calculation (Eq. 10).
SET_SPOT_AREA	Set the filter spot area (Eq. 6) in m ² .
SET_TIME_ZONE	Set time zone
SPD	Set the pump duty cycle to change the instrument flow.

5.3. Action command

Command	Description
CHANGE_FILTER	<i>Open the measuring chamber, filter advancement and close the measuring chamber. Command can only be run when the measurement is not active, after this action measure does not start automatically.</i>
CLOSE_HEAD	<i>Close measurement chamber.</i>
DELETE_ALL_MEASURES	<i>Delete all the measures stored in the SD card.</i>
DELETE_DATE_MEASURES	<i>Delete the measures of the given day stored in the SD card.</i>
DELETE_LOGS	<i>Remove all system logs.</i>
DISABLE_FLUXMETER	<i>Disable the reading of the fluxmeter, to use only if the fluxmeter is broken and an external fluxmeter is used. Data are given with a default flux and must be corrected during post-analysis.</i>
DISABLE_PUMP	<i>This command allows to perform measures without flow (pump off). Data are then given with a default flow . Used mainly for electronic noise check. Command can only be run when the measurement is not active, after this action measure does not start automatically.</i>
DISABLE_REMOTE_CONSOLE	<i>Disable the telnet communication.</i>
ENABLE_FLUXMETER	<i>Enable the reading of the fluxmeter (default condition).</i>
ENABLE_PUMP	<i>Enable pump, to use after the disabled pump command to return in normal operation mode. Command can only be run when the measurement is not active, after this action measure does not start automatically.</i>
ENABLE_REMOTE_CONSOLE	<i>Enable the telnet communication.</i>
FORWARD_REEL	<i>Advancing the filter. This command can be executed only when the measurement chamber is open.</i>
LED_BLUE	<i>Activate only blue led. To be used in calibration mode.</i>
LED_GREEN	<i>Activate only green led. To be used in calibration mode.</i>
LED_IR	<i>Activate only ir led. To be used in calibration mode.</i>
LED_NONE	<i>None of the LED is active. To be used in calibration mode for dark current.</i>
LED_RED	<i>Activate only red led. To be used in calibration mode.</i>
LED_UV	<i>Activate only uv led. To be used in calibration mode.</i>
OPEN_HEAD	<i>Open the measurement chamber. Command can only be run when the measurement is not active.</i>
START_CALIBRATION	<i>Put the instrument in calibration mode. Command can only be run when the measurement is not active.</i>
START_MEASURE	<i>Start measures.</i>
START_PUMP	<i>Activate the pump.</i>
STOP_CALIBRATION	<i>Stop the calibration mode.</i>
STOP_MEASURE	<i>Stop the measure</i>
STOP_PUMP	<i>Stop the pump</i>

5.4. Other commands

Command	Description
ADMIN	<i>Enable the administration mode. Reserved for the producer – need a password.</i>
EXIT	Only over TCP/IP. <i>Disable remote console (telnet).</i>

6. Changing the tape filter

1. Stop measures
2. Open head
3. Open the Plexiglas cover
4. Advance the remaining filter by manually running the right knob
5. When the filter is completely removed, unscrew the right knob and remove the used tape filter from the support
6. Place the right knob back on the support
7. Unscrew the left knob and insert the new filter tape on the support
8. Block the filter screwing back the knob, insert the filter tape from the left side of the control camera and pass it through, then stop it. Use the adhesive tape on the right support
9. Close the Plexiglas cover
10. Close head
11. Advance the filter (forward_reel) to check if the operation was successful

Figure 3 shows the correct position of the tape filter inside the instrument.

APPENDICE A: data format

In the following table is reported the data format obtained in streaming or by using GET_DATE_MEASURES, GET_DATE_RANGE_MEASURES and GET_LAST_MEASURES commands.

<i>Parameter</i>	<i>header</i>	<i>Unit</i>	<i>Parameter</i>	<i>header</i>	<i>Unit</i>
Date	DATE	yyyy-mm-dd	UV transmittance	TRANS_UV	
Time	TIME	HH:MM:SS	IR absorption coefficient	KABS_IR	m ⁻¹
Decimal day (0 = 1 st January)	NUMERIC_DAY	days	RED absorption coefficient	KABS_RED	m ⁻¹
Venturi tube differential pressure	FLUX_PRESS	mbar	GREEN absorption coefficient	KABS_GREEN	m ⁻¹
Flow	FLUX_L/M	l/min	BLUE absorption coefficient	KABS_BLUE	m ⁻¹
LED temperature	LED_TEMP	°C	UV absorption coefficient	KABS_UV	m ⁻¹
Air temperature (in Venturi tube)	AIR_TEMP	°C	Equivalent Black carbon IR	BC_USER	µg/m ³
Main-Board temperature	INT_TEMP	°C	Ambient pressure	ENVIRO_PRESS	hPa
IR transmittance	TRANS_IR		Flags	FLAGS	hex
RED transmittance	TRANS_RED		AAE 870-420	AAE	
GREEN transmittance	TRANS_GREEN		RH	AIR_HUMI	%
BLUE transmittance	TRANS_BLUE		Time Zone	TIME_ZONE	
			SMA	SMA	minutes



APPENDIX B: Flags

The last field in the data (Flags) is an hexadecimal number that report system status.
The table below shows the value of the flags and their meaning.

<i>Flag</i>	<i>Meaning</i>
1	Cooler fan active
2	Pump disabled.
4	Start of new measure

In case more flag are detected together they are added in the flags. For example, flag 3 is the sum of flag 1 and 2, that means that the cooler fan is active and there is a pump disabled.

APPENDIX C: Electrical and mechanical specification

Electrical		
<i>Power supply</i>	<i>220 Vac</i>	<i>Max 0.5A</i>
<i>Power consumption</i>		<i>36 Watt</i>
<i>Mains filter</i>	<i>Arctronics AR3 series</i>	
<i>Protection</i>	<i>2 x fuse (5x20)</i>	<i>2A</i>
<i>Power plug</i>	<i>IEC</i>	
<i>Ethernet</i>	<i>RJ45</i>	<i>10-100 Mbps</i>
<i>RS232</i>	<i>DB9 (3 pin)</i>	<i>57.600 Kbps</i>

Mechanical		
<i>Dimension</i>	<i>19" x 6 Unit</i>	<i>Max 460 x 300 x 270 mm</i>
<i>Weight</i>	<i>11 Kg</i>	
<i>Tape filter</i>	<i>width 40 mm</i>	<i>type GF10</i>
<i>Air Inlet</i>	<i>D= 6 mm</i>	
<i>Air outlet</i>	<i>D= 6 mm</i>	
<i>Air flow (adjustable)</i>	<i>1 – 2.5 l/min</i>	
<i>Working temperature</i>	<i>10-30 C°</i>	



DBAP5

Manual version 1.16-EN
Firmware 1.5.2 or higher

APPENDIX D: Normative

The instrument is harmonized with the following standards:

- CE, UL60950-1, EN60950-1
- CEI EN 61010-1
- 2006/95/EC
- 2004/108/CE
- 1990-03-05/4690 /art 3 22012008/37

Intended use: Laboratory equipment for environmental monitoring.

Via S. Andrea di Rovereto 33A CS - 16043 Chiavari (GE) Italy - Tel. +39 0185 318444

Email: len@len.it url: <http://www.len.it>

APPENDIX E: Poster PM2022

Development of a new photometer for measuring the absorption coefficient at 5 different wave length Dual Beam Absorption Photometer 5 (DBAP5), description and first comparisons

MAURIZIO BUSETTO¹, LAURA RENZI¹, LUCA DI LIBERTO² AND ANGELO LUPI³
¹ Istituto di Scienze dell'Atmosfera e del Clima, Consiglio Nazionale delle Ricerche, Bologna
² Istituto di Scienze dell'Atmosfera e del Clima, Consiglio Nazionale delle Ricerche, Roma
³ Istituto di Scienze Polari, Consiglio Nazionale delle Ricerche, Bologna




Motivation DBAP5 description Instrumental intercomparison

Black Carbon (BC) is the main absorbing component of atmospheric aerosol. Generated by the incomplete combustion of fuels, biomass burning, internal combustion engines (especially diesel), and industrial processes, it has a lifetime in the atmosphere of days to weeks.

By strongly absorbing solar radiation and reducing albedo, when deposited on snow and ice, BC has a warming effect on climate [1], the quantification of which is still uncertain.


Moreover, epidemiological studies provide evidence of the association of cardiopulmonary morbidity and mortality with exposure to black carbon.

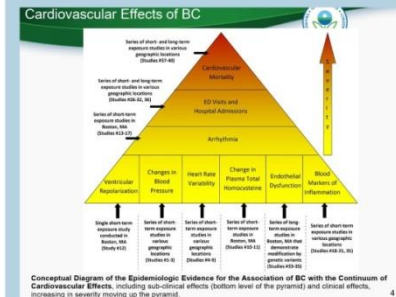

These features have made BC an important pollutant to monitor, understand, and manage despite not currently having direct legal limits imposed in Europe and elsewhere (Reche et al., 2011; European Environment Agency, 2016).

Absorption coefficient Spectra Dependence (AAE)


Black carbon is not the only absorbing component of atmospheric aerosol, dust and organic compounds also absorb the radiation but with a different spectral dependence of the absorption coefficient, quantified by the Absorption Angstrom Exponent (AAE).

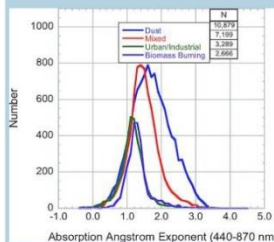


This is thus a key parameter to evaluate aerosol composition and sources.

D. M. Gilles et al. 2012

The most popular method for measuring BC concentration and spectrally resolved aerosol absorption coefficient is the filter-based absorption photometer. Particles are collected on filter and as the particle load increase, the light passing through the filter decrease. DBAP5 is a 5 wavelength (from 420 to 870 nm, filter-based photometer based on the dual beam technology (Dual Beam Absorption Photometer).

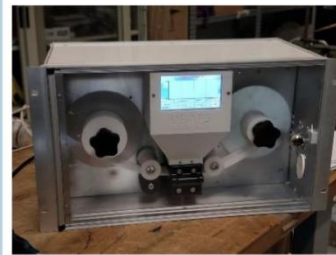

Measuring the variation in filter light transmission at increasing particle load it evaluate the attenuation coefficients, then converted in the absorption coefficient by applying the appropriate filter correction equations.

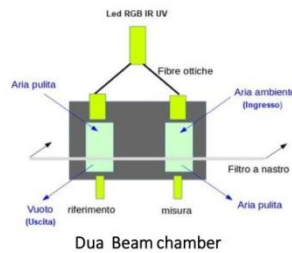


Equivalent Black Carbon concentration can be then derived using the appropriate Mass Absorption Coefficients. The dual beam technology is a technique that simultaneously compare the absorption of the particulate matter with the absorption on the white filter, providing more precise measurement especially in low concentration site.

Technica Specification

LED	870,634,522,465,420 nm
Air flow	1-5 l/m
Filter tape	Glass fiber GF-10 w=40mm
communication	RS232, telnet

Urban site Bo gna with an AE33 (spring 2021)


The BC [ng/cm³] DBAP5 has been evaluated from absorption coefficient at 870 nm

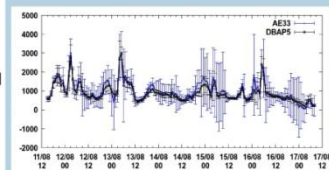

One hour averaged data has been used to compare DABP5 and AE33 measures. Correlation is very good, Comparing the SD, the


DBAP5's SD is generally lower than AE33's, indicating less noise in the DBAP5 measurements in Urban areas

During this period AAE has been evaluated from DBAP5 data, results agree with [3], showing typical values of urban sites

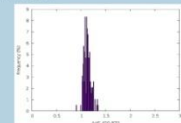

Rura site San Pietro Capofiume, with a MAAP, Spring 2022

The BC for DBAP5 has been evaluated comparing it with BC measured using a Multi Angle Absorption Photometer.

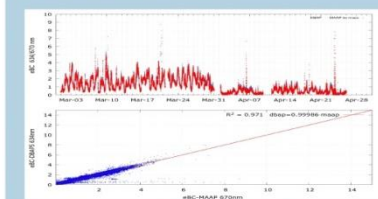

One-minute data have been compared for two months. The results show a powerful agreement between two instrumental setup.

- 1) Reche, C. et al.: New considerations for PM, Black Carbon and Atmos. Meas. Tech.
- 2) European Environment Agency: Black carbon: Better monitoring needed to assess health and climate change impacts
- 3) An analysis of AERONET aerosol absorption properties and classifications representative of aerosol source regions. D. M. Gilles et al. 2012. J. Geophys. Res.
- 4) S. K. Grange et al.: Evaluation of EBC source apportionment using observations from Switzerland 1885 particle number concentration for air quality monitoring across different European cities